




Scan-RAM 2™

Calibration Sources


www.lablogic.com

 **LabLogic**
EXPERIENCE & EXPERTISE


Scan-RAM 2 - Three source calibration bed

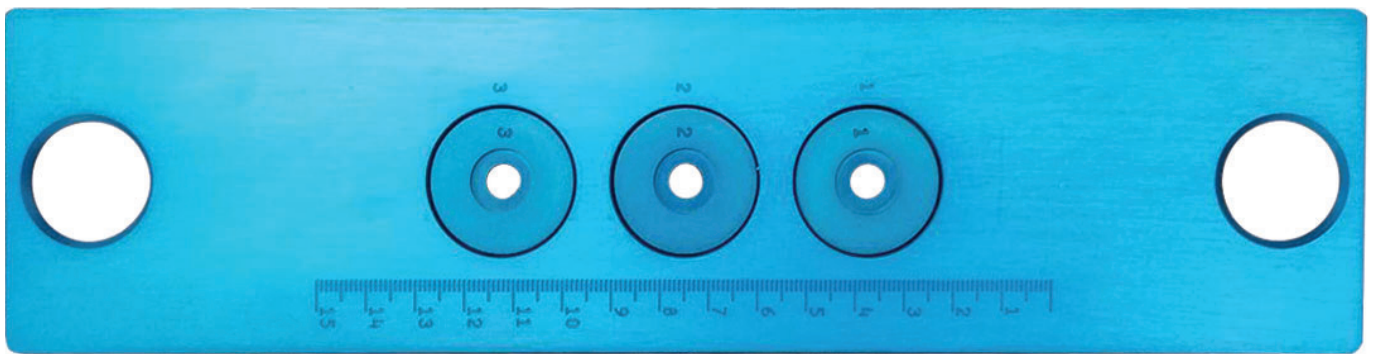

Scan-RAM 2 - Four source calibration bed

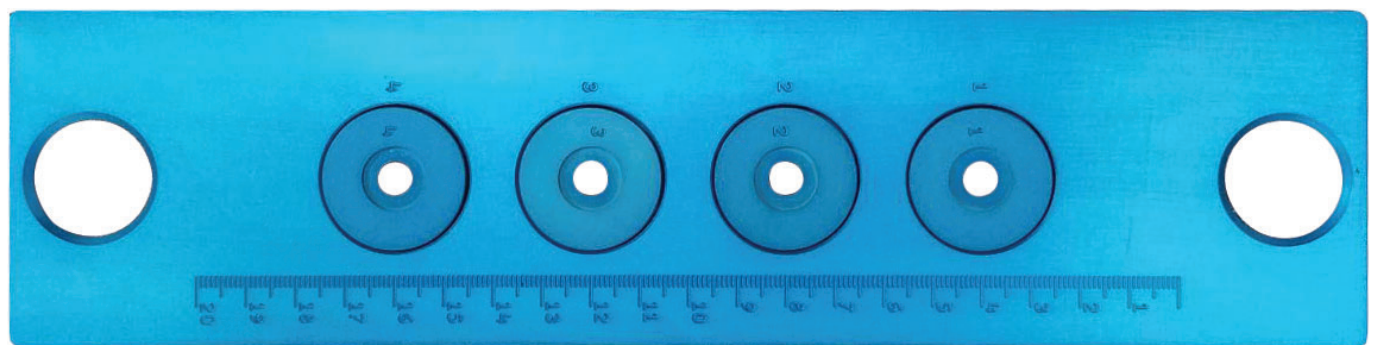


NaI SiPM with Co-57

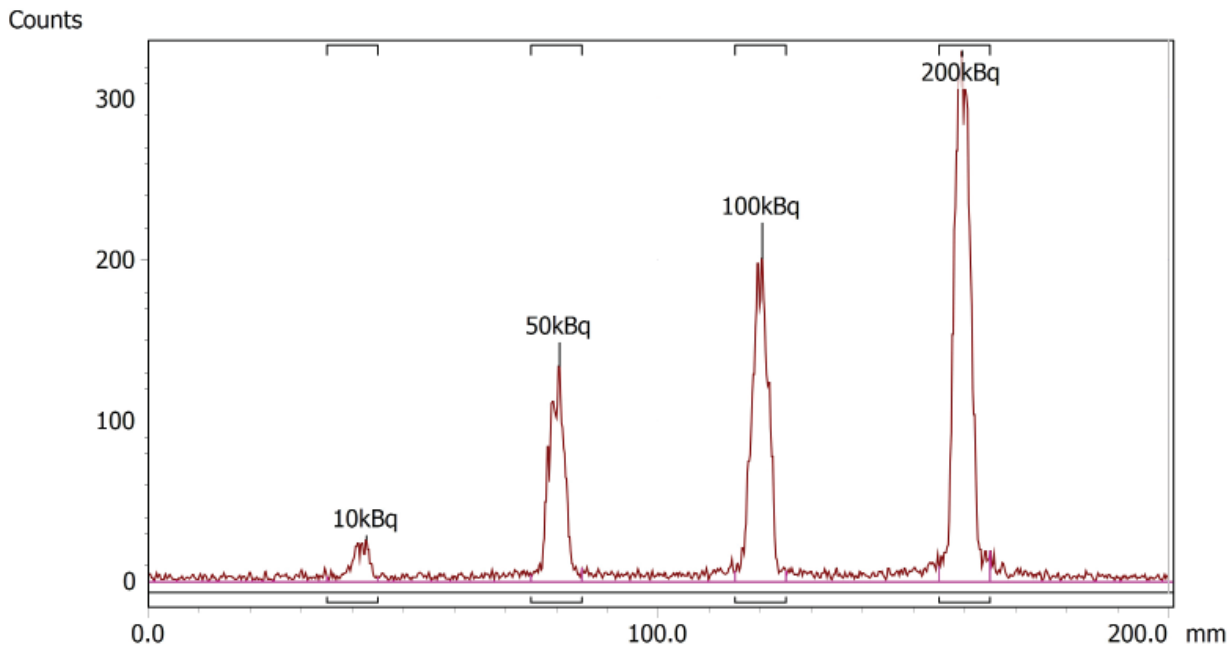


Name	Start (mm)	End (mm)	Retention (RF)	Area (Counts)	% ROI (%)	% Total (%)
10kBq	35.0	45.0	0.214	466	3.86	3.01
50kBq	75.0	85.0	0.403	2207	18.27	14.27
100kBq	115.0	125.0	0.602	3575	29.59	23.12
200kBq	155.0	165.0	0.798	5832	48.28	37.71
4 Peaks				12080	100.00	78.11

NaI SiPM with Cs-137

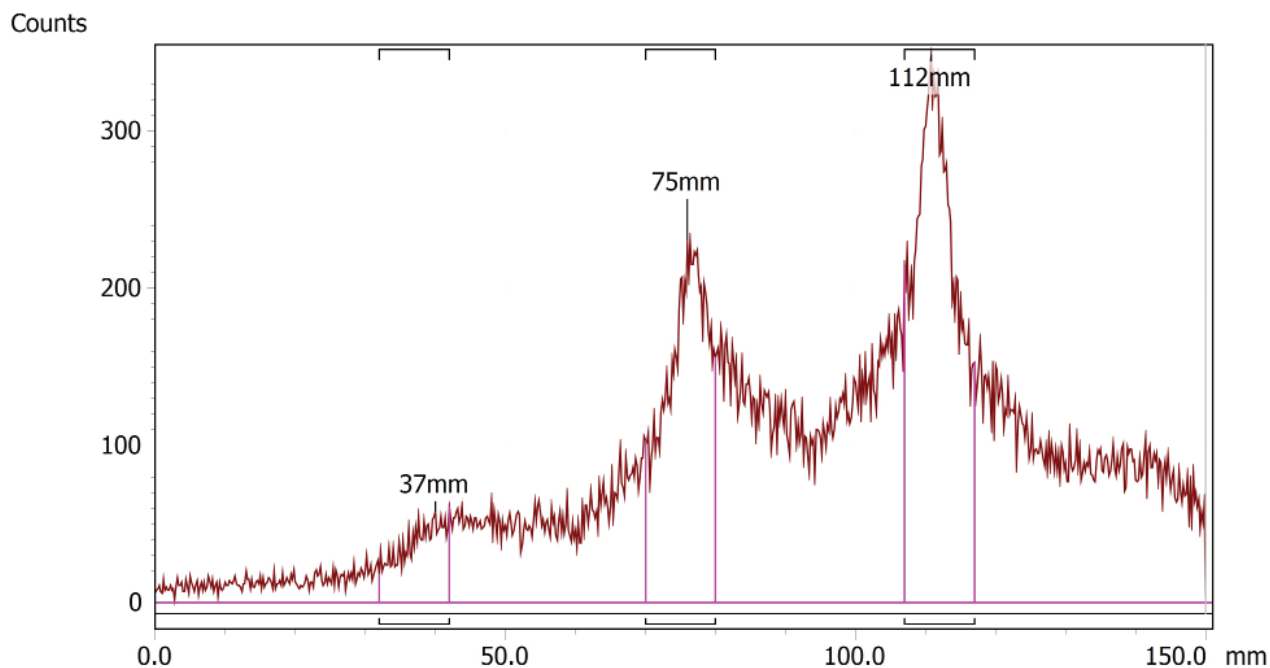


Name	Start (mm)	End (mm)	Retention (RF)	Area (Counts)	% ROI (%)	% Total (%)
37mm	32.0	42.0	0.267	1899	8.67	2.90
75mm	70.0	80.0	0.507	8039	36.72	12.28
112mm	107.0	117.0	0.739	11956	54.61	18.26
3 Peaks				21894	100.00	33.44

Plastic SiPM with Co-57

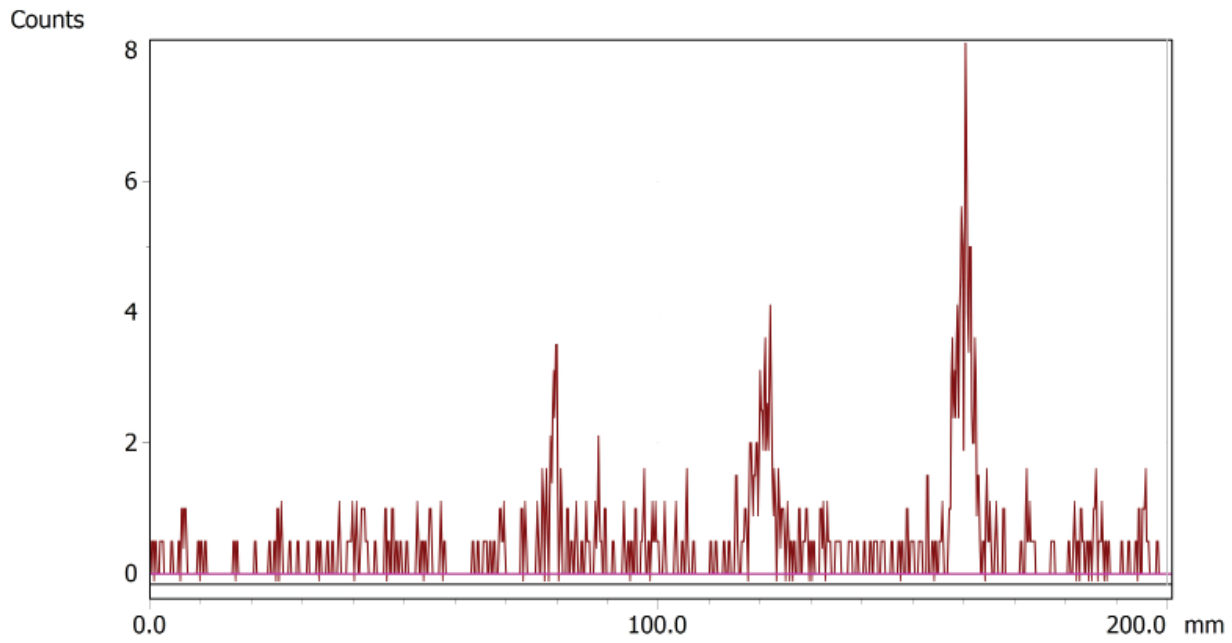

Name	Start (mm)	End (mm)	Retention (RF)	Area (Counts)	% ROI (%)	% Total (%)
0 Peaks				0	100.00	0.00

Plastic SiPM with Cs-137

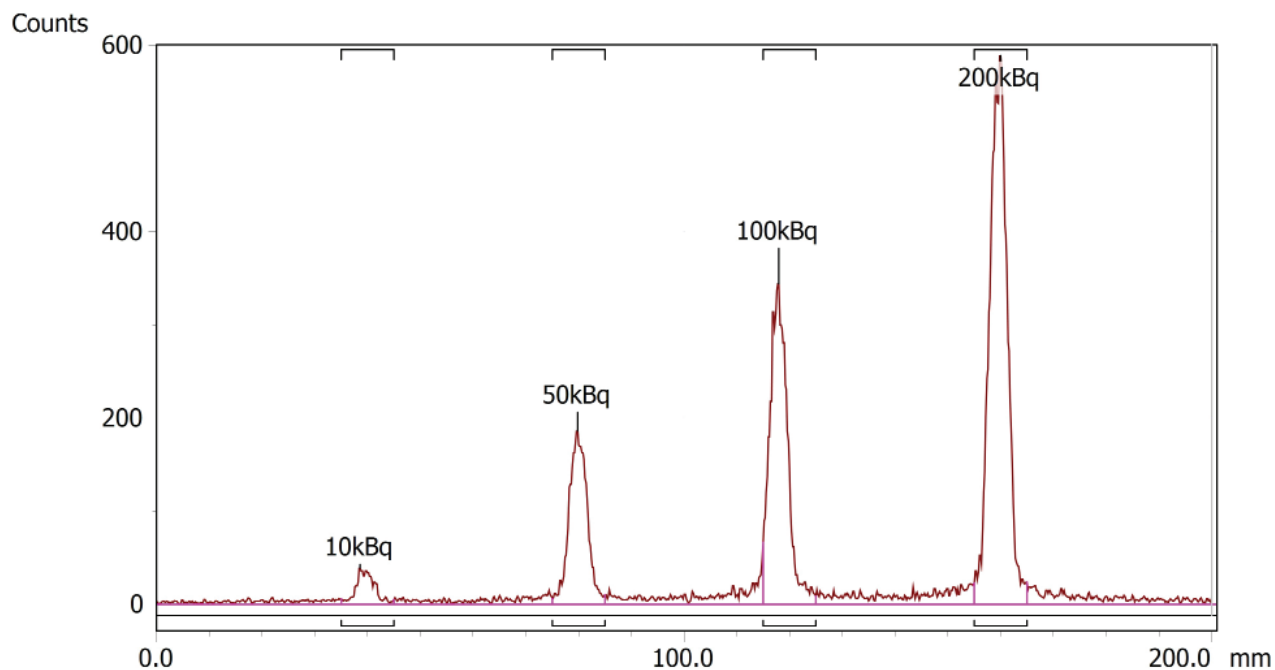

Name	Start (mm)	End (mm)	Retention (RF)	Area (Counts)	% ROI (%)	% Total (%)
10kBq	35.0	45.0	0.193	703	3.27	2.60
50kBq	75.0	85.0	0.399	3489	16.23	12.92
100kBq	115.0	125.0	0.590	6324	29.42	23.41
200kBq	155.0	165.0	0.800	10981	51.08	40.65
4 Peaks				21497	100.00	79.57

Service and Support

Users of our systems can benefit from our comprehensive, fully inclusive service and support.

We can give reassurance that if things go wrong or you need expert advice, help is only an e-mail or phone call away.



Validation Services

Our Validation Service enables you to implement and get maximum value from your investments as soon as possible.

We work as a partner with your Quality Manager, System Manager and users to provide a tailored Validation Plan, suited to your needs. Our Validation Specialists who have many years' of experience in GLP system validation, detailed knowledge of our systems, together with other industry standard systems to help you meet company and regulatory requirements.

Training

LabLogic can provide a variety of training courses and workshops to help you get the most out of your instrument and software.

All training is performed by our expert Product and Support Specialists who have many years experience in the development and use of the instruments and software.

Certificates can be provided to complement your internal GLP training records.

Europe & Worldwide

LabLogic Systems Limited

Innovation House, 6 Europa View
Sheffield, S9 1XH, UK

E-mail: solutions@lablogic.com

Tel: +44 (0)114 266 7267

www.lablogic.com

USA & Canada

LabLogic Systems, Inc.

3901 Centerview Drive, Suite B
Chantilly, VA 20151, USA

E-mail: solutions@lablogic.com

Tel: +1-703-429-4209

www.lablogic.com



Certificate No: 1535
ISO 9001



Certificate No: 10926
ISO 9001