

Marc 7000

Tritium Air Sampler



The Marc 7000 tritium sampler from SDEC is the perfect instrument for measuring low levels of tritium in air.

- Trapping yield of 99% by bubbling air through demineralised water.
- Gas (HT) and vapor (HTO) tritium forms can both be collected with the catalytic oven.
- Reduced evaporation due to the cooled trapping system, allowing weekly collection.
- Easy-to-use, with instant opening cabinet for sample retrieval.
- Accurate – electronic, accredited airflow meter.

 **LabLogic**

EXPERIENCE & EXPERTISE

Marc 7000

The Marc 7000 is widely used and recognised within the international nuclear industry, and in particular; nuclear power plants, nuclear research centres, radioactive waste treatment facilities and isotope laboratories.

The bubbler has been specifically designed with efficient tritium capture in mind, using a series of four vials, a cooling system and a catalytic oven to collect both tritium vapor (HTO) and gaseous tritium (HT).

The tritium activity in the collected sample can be measured with a liquid scintillation counter on a daily, weekly or monthly basis, and can then be used in combination with the sampled air volume to calculate the tritium-in-air concentration. This provides an efficient method of monitoring tritium levels with a much higher sensitivity than even the most sophisticated real-time tritium monitor.

Airflow process

A pre-filter paper in the inlet (Ø 45 mm) prevents dust intake and the electronic flow meter is protected by Gortex® filters. The airflow passes into glass vials (250 ml capacity) through stainless steel air tubing. The air flow can be set from 10 to 50 litres per hour, regulated by a certified airflow meter.

Oxidation oven

Stainless steel tubular oven equipped with Palladium alumina catalyst pellets. The oven temperature can be set between +200°C and +500°C.

Cooling system of the collection vials

The sampler is fitted with a condenser cooling block, which allows the vials to be cooled to between +5°C and +15°C (depending on the ambient temperature). A pump ensures flow of the cooling liquid and a level gauge allows the direct control of the liquid level in the circuit. All tubing is made of stainless steel.

Alarms and defaults

- Alarm buzzer for all detected defaults.
- Memory and recovery of the last 8 defaults via RS-232.

Efficiency

HTO trapping yield	96% ± 4%
Oven conversion yield	95% ± 4%
Reproducibility of the airflow	± 0.8%
Airflow accuracy	± 1%

No response to other radioactive elements.

Options

- Condensation collector tray (recommended).
- Alarm state relay.
- J-bus protocol for remote control.
- External presurre regulator.
- Flashing light alarm signal.

Specification

Display	LCD display featuring: Oven temperature, cooling liquid temperature. Instant air flow and total volume. Duration of sampling and alarm.
Front Panel	Scratch proof lexan cover.
Frame	Monocoque in aluminium alloy. Decontamination compliant housing plant
Power Supply	230 V / 50Hz or 120 V / 60 Hz IEC plug.
Power	700 Watts max.
Regulation	Electronic control of the airflow, oven temperature, cool temperature.
Electrical Protection	Differential circuit breaker (sensitivity = 30 mA).
Inlet Connections	Flexible plastic tube inner diameter 6 mm with fastenings.
Temp. (Operating)	+ 2°C to + 45°C
Temp. (Storage)	-5°C to + 70°C
Dimensions	27.5" x 10" x 10.5" (700 x 265 x 270 mm)
Required Space	39" x 23.5" x 20.5" (1000 x 600 x 530 mm)
Weight	64 lbs (29 kg)

USA & Canada

LabLogic Systems, Inc.

East Pointe Park, 1040 East Brandon Blvd.
Brandon, FL 33511-5509, USA

E-mail: solutions@lablogic.com

Tel: +1-813-626-6848

Fax: +1-813-620-3708

Web: www.lablogic.com



Certificate No: 10926
ISO 9001

Europe & Worldwide

LabLogic Systems Limited

Paradigm House, 3 Melbourne Avenue
Broomhill, Sheffield, S10 2QJ, UK

E-mail: solutions@lablogic.com

Tel: +44 (0)114 266 7267

Fax: +44 (0)114 266 3944

Web: www.lablogic.com



Certificate No: 1535
ISO 9001



EXPERIENCE & EXPERTISE